


Preparation of CRM samples in powder format

1

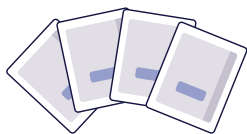

Check the certificate to select desired level of microorganisms.

2



Accurately weigh a minimum of 10 g of powder.

Prepare your desired dilution, any ratio can be used, e.g. 1:10.

The material is now ready for testing.

3

There are several options on how to use the resuscitated material

A



(Dilute and) Directly plate onto agar/petrifilm or other.

B



Weigh and dilute a new sample matrix of your choice. Add desired volume from your primary diluted CRM (from step 2).

C



Prepare a 1 liter water sample. Add desired volume from your primary diluted CRM (from step 2).

4



Proceed to testing according to your method of choice.

More information



eurofins.no/CRM